

Now please try to solve the problem according to the instructions below.

Keep pressing MODE key and it will display the using segments of the display screen after about 3 seconds.

Unloos the MODE key and then press CAL key it will display line 0.

Now confirm that there is nothing on the pan.

Press TARE key the balance system will start to collect the data of the zero point.

After then I'll prompt the calibration weight which should be put on the pan.

Lin 1 1kg

Lin 2 2kg

Lin 3 3kg

Lin 4 4kg

Lin 5 5kg

After putting the corresponding weight press TARE key for the confirmation.

After then, the system will collect all the corresponding data and it will display the final group of the weighing data.

After the linear adjustment you can check the linear error.

If it's not eligible you can adjust the linear again.

Avoid the error influence of the artificial operation.

That's all.